

# Product Sheet

## H\_CD25 CD122 CD132 Reporter Cell Line


Catalog number: GM-C29055

Version 3.3.1.241219

Interleukin-2 (IL-2) is an important cytokine that mainly plays a regulatory role in the immune system. It works by binding to IL-2 receptors on the surface of lymphocytes, activating a series of signaling pathways that influence the proliferation and activity of immune cells.

The IL-2 receptor has two forms: a heterotrimer composed of IL-2R $\alpha$ , IL-2R $\beta$ , and IL-2R $\gamma$ , and a heterodimer composed of IL-2R $\beta$  and IL-2R $\gamma$ . When IL-2 binds to its receptor, it activates the JAK signaling pathway, which in turn activates the transcription factor STAT. The phosphorylated STATs form dimers or tetramers and move into the cell nucleus, regulating the expression of specific genes to promote the immune response.

H\_CD25 CD122 CD132 Reporter Cell Line is a clonal stable cell line constructed using lentiviral technology, constitutive expression of the IL-2R $\alpha$ (CD25), IL-2R $\beta$ (CD122) and IL-2R $\gamma$ (CD132) gene, along with signal-dependent expression of a luciferase reporter gene. When IL-2 binds to IL-2R, it activates downstream signaling pathways, leading to the expression of luciferase. The luciferase activity measurement indicates the activation level of the signaling pathway and can thus be used to evaluate the in vitro effects of drugs related to IL-2.

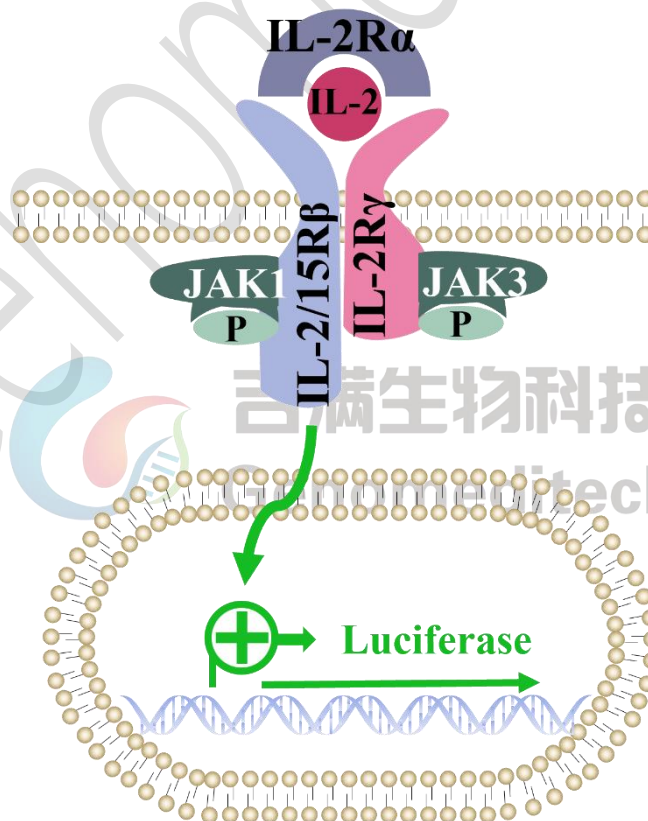

## Specifications

<b>Quantity</b>	5E6 Cells per vial, 1 mL
<b>Product Format</b>	1 vial of frozen cells
<b>Shipping</b>	Shipped on dry ice
<b>Storage Conditions</b>	Liquid nitrogen immediately upon receipt
<b>Recovery Medium</b>	RPMI 1640+10% FBS+1% P.S
<b>Growth medium</b>	RPMI 1640+10% FBS+1% P.S+3.5 µg/mL Blasticidin+400 µg/mL G418+200 µg/mL Hygromycin+0.75 µg/mL Puromycin
<b>Note</b>	None
<b>Freezing Medium</b>	90% FBS+10% DMSO
<b>Growth properties</b>	Suspension
<b>Growth Conditions</b>	37°C, 5% CO <sub>2</sub>
<b>Mycoplasma Testing</b>	The cell line has been screened to confirm the absence of Mycoplasma species.
<b>Safety considerations</b>	Biosafety Level 2
<b>Note</b>	It is recommended to expand the cell culture and store a minimum of 10 vials at an early passage for potential future use.

## Materials

Reagent	Manufacturer/Catalogue No.
RPMI 1640	VivaCell/C3010-0500
Fetal Bovine Serum	Cegrogen biotech/A0500-3010
Pen/Strep	Thermo/15140-122
Blasticidin	Genomeditech/ <a href="#">GM-040404</a>
G418	Genomeditech/ <a href="#">GM-040402</a>
Hygromycin	Genomeditech/ <a href="#">GM-040403</a>
Puromycin	Genomeditech/ <a href="#">GM-040401</a>
H_CD122 CD132 Reporter Cell Line	Genomeditech/ <a href="#">GM-C29054</a>
Recombinant Human IL-2	Novoprotein/C013
Anti-CD25 hIgG1 Antibody(Basiliximab)	Genomeditech/ <a href="#">GM-52329AB</a>
Anti-CD122 hIgG1 Antibody(HuABC-2)	Genomeditech/ <a href="#">GM-52319AB</a>
Anti-CD132(IL2RG) hIgG4 Antibody(REGN7257)	Genomeditech/ <a href="#">GM-52334AB</a>
GMOne-Step Luciferase Reporter Gene Assay Kit	Genomeditech/ <a href="#">GM-040503</a>

## Figures

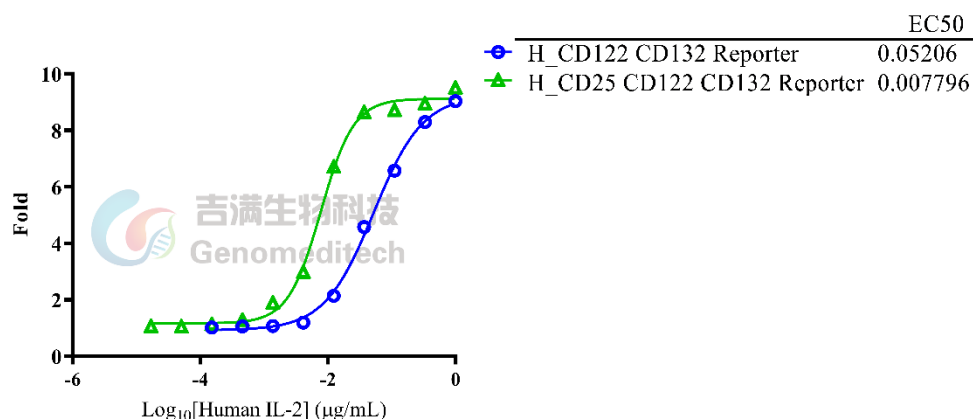


Figure 1 | Response to Recombinant Human IL-2. The H\_CD122 CD132 Reporter Cell Line (Cat. [GM-C29054](#)) and H\_CD25 CD122 CD132 Reporter Cell Line (Cat. [GM-C29055](#)) at a concentration of 1E5 cells/well (96-well format) were stimulated with serial dilutions of Recombinant Human IL-2 (Novoprotein/C013) in assay buffer (RPMI 1640 + 1% FBS + 1%P.S) for 16 hours. The firefly luciferase activity was measured using the GMOne-Step Luciferase Reporter Gene Assay Kit (Cat. [GM-040503](#)). The maximum induction fold was approximately [9.0] and [9.5], respectively. Data are shown by drug mass concentration.

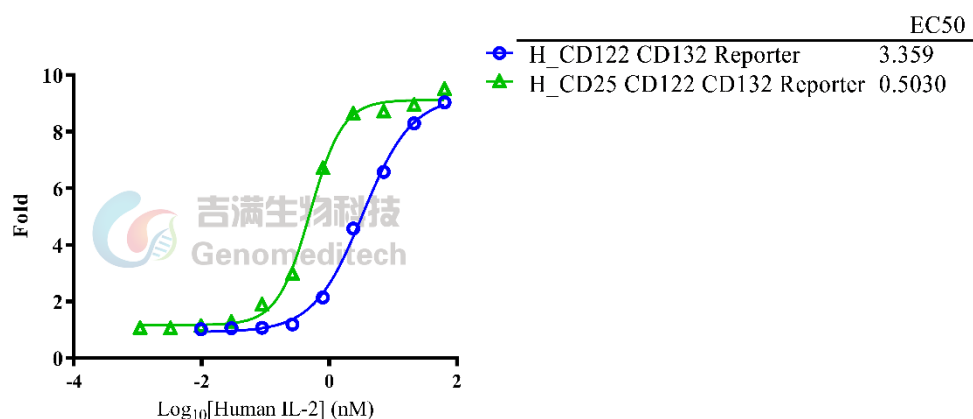


Figure 2 | Response to Recombinant Human IL-2. The H\_CD122 CD132 Reporter Cell Line (Cat. [GM-C29054](#)) and H\_CD25 CD122 CD132 Reporter Cell Line (Cat. [GM-C29055](#)) at a concentration of 1E5 cells/well (96-well format) were stimulated with serial dilutions of Recombinant Human IL-2 (Novoprotein/C013) in assay buffer (RPMI 1640 + 1% FBS + 1%P.S) for 16 hours. The firefly luciferase activity was measured using the GMOne-Step Luciferase Reporter Gene Assay Kit (Cat. [GM-040503](#)). The maximum induction fold was approximately [9.0] and [9.5], respectively. Data are shown by drug molar concentration.

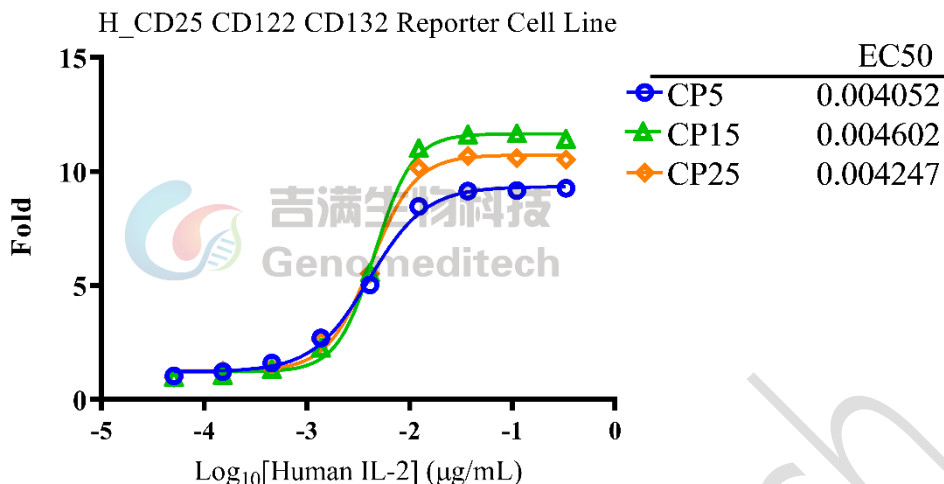


Figure 3 | The passage stability of response to Recombinant Human IL-2. The passage 5, 15 and 25 of H\_CD25 CD122 CD132 Reporter Cell Line (Cat. GM-C29055) at a concentration of 1E5 cells/well (96-well format) was stimulated with serial dilutions of Recombinant Human IL-2 (Novoprotein/C013) in assay buffer (RPMI 1640 + 1% FBS + 1%P.S) for 16 hours. The firefly luciferase activity was measured using the GMOne-Step Luciferase Reporter Gene Assay Kit (Cat. GM-040503). Data are shown by drug mass concentration.

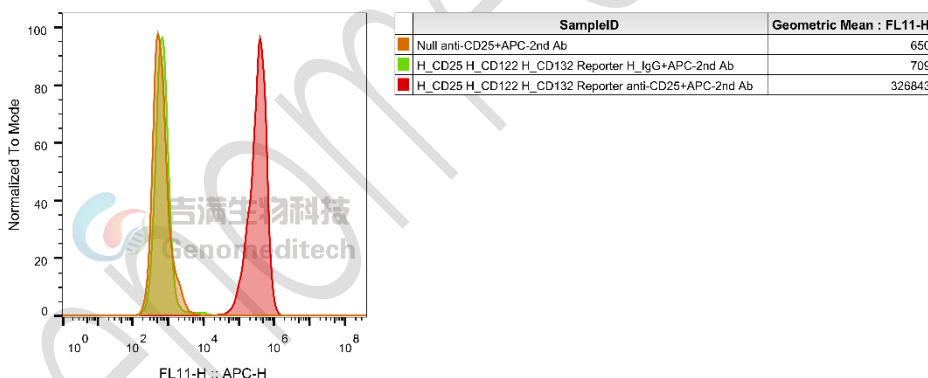


Figure 4 | H\_CD25 CD122 CD132 Reporter Cell Line (Cat. GM-C29055) was determined by flow cytometry using Anti-CD25 hIgG1 Antibody(Basiliximab) (Cat. GM-52329AB).

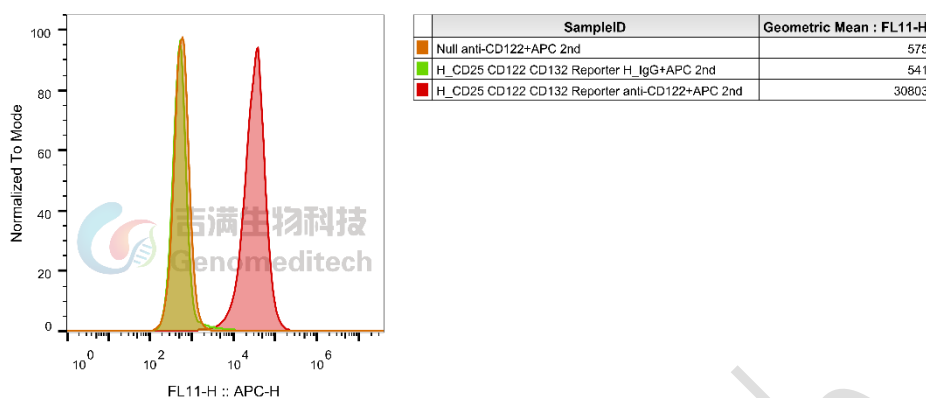


Figure 5 | H\_CD25 CD122 CD132 Reporter Cell Line (Cat. GM-C29055) was determined by flow cytometry using Anti-CD122 hlgG1 Antibody(HuABC-2) (Cat. [GM-52319AB](#)).

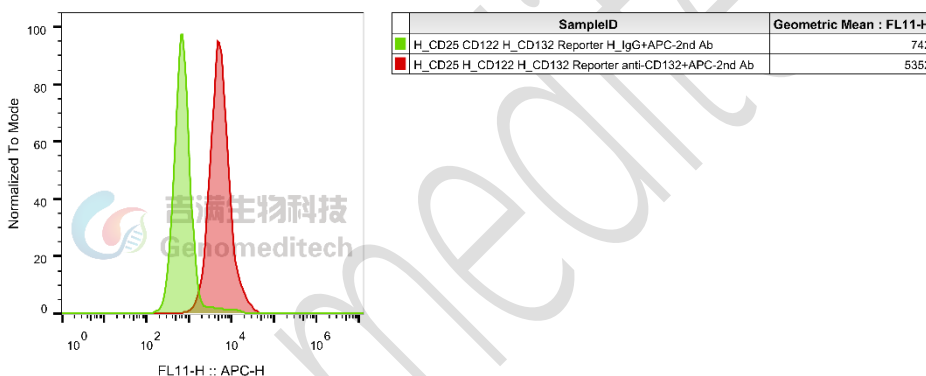


Figure 6 | H\_CD25 CD122 CD132 Reporter Cell Line (Cat. GM-C29055) was determined by flow cytometry using Anti-CD132(IL2RG) hlgG4 Antibody(REGN7257) (Cat. [GM-52334AB](#)).

## Cell Recovery

Recovery Medium: RPMI 1640+10% FBS+1% P.S

To insure the highest level of viability, thaw the vial and initiate the culture as soon as possible upon receipt. If upon arrival, continued storage of the frozen culture is necessary, it should be stored in liquid nitrogen vapor phase and not at -70°C. Storage at -70°C will result in loss of viability.

- Thaw the vial by gentle agitation in a 37°C water bath. To reduce the possibility of contamination, keep the O-ring and cap out of the water. Thawing should be rapid (approximately 2 - 3 minutes).
- Remove the vial from the water bath as soon as the contents are thawed, and decontaminate by dipping in or spraying with 70% ethanol. All of the operations from this point on should be carried out under strict aseptic conditions.
- Transfer the vial contents to a centrifuge tube containing 5.0 mL complete culture medium. And spin at approximately 176 x g for 5 minutes. Discard supernatant.

- d) Resuspend cell pellet with the recommended complete medium. And dispense the suspension into 1 - 2 T-25 culture flasks.
- e) Incubate the culture at 37°C in a suitable incubator. A 5% CO<sub>2</sub> in air atmosphere is recommended if using the medium described on this product sheet.

## Cell Freezing

Freezing Medium: 90% FBS+10% DMSO

- a) Centrifuge at 176 x g for 3 minutes to collect cells.
- b) Resuspend the cells in pre-cooled freezing medium and adjust the cell density to 5E6 cells/mL.
- c) Aliquot 1 mL into each vial.
- d) Place the vial in a controlled-rate freezing container and store at -80°C for at least 1 day, then transfer to liquid nitrogen as soon as possible.

## Cell passage

Growth medium: RPMI 1640+10% FBS+1% P.S+3.5 µg/mL Blasticidin+400 µg/mL G418+200 µg/mL Hygromycin+0.75 µg/mL Puromycin

Approximately 48-72 hours after the initial thawing, the cells can be passaged for the first time. After this initial passage, the culture medium can be adjusted to growth medium supplemented with antibiotics. If cells are not passaged within 48 hours, it is recommended to add some fresh recovery medium and place the flask horizontally.

- a) When the cell density reaches 1.5 - 2E6 cells/mL, subculture the cells. Do not allow the cell density to exceed 2E6 cells/mL.
- b) It is recommended to use T-25 flasks for subculturing.
- c) These cells are suspension cells, and it is recommended to use the "half-medium change" method to maintain optimal cell conditions during passaging.
- d) During passaging, you can directly add fresh growth medium to the culture flask, gently pipette to resuspend the cells, and then transfer the cell suspension to a new T-25 flask for continued culture.

**Subcultivation Ratio: Maintain cultures at a cell concentration between 3E5 and 1E6 viable cells/mL.**

**Medium Renewal: Every 2 to 3 days**

## Notes

- a) These cells are sensitive to density, so please ensure that the cell density is maintained within an appropriate range during culture and subculturing.
- b) During the first passage, pay attention to the nutrient supply; if not subculturing, make sure to add fresh recovery medium every other day as needed.

## Related Products

IL-15

H_IL15 Reporter Cell Line	Cynomolgus_CD122 HEK-293 Cell Line
H_CD122 CHO-K1 Cell Line	H_CD122 HEK-293 Cell Line
H_CD215(IL15RA) HEK-293 Cell Line	
IL-2	
H_CD122 CD132 Reporter Cell Line	H_IL2 Reporter Cell Line
H_IL2 Reporter DDX35TM Cell Line	Cynomolgus_CD25 HEK-293 Cell Line
H_CD25 CHO-K1 Cell Line	H_CD25 HEK-293 Cell Line
Anti-CD122 hIgG1 Antibody(HuABC-2)	Anti-CD132(IL2RG) hIgG4 Antibody(REGN7257)
Anti-CD25 hIgG1 Antibody(Basiliximab)	Anti-mouse CD25 mIgG2a Antibody(PC-61.5.3)
Anti-mouse CD25 RIgG1 Antibody(PC-61.5.3)	

## Limited Use License Agreement

Genomeditech (Shanghai) Co., Ltd grants to the Licensee all intellectual property rights, exclusive, non-transferable, and non-sublicensable rights of the Licensed Materials; Genomeditech (Shanghai) Co., Ltd will retain ownership of the Licensed Materials, cell line history packages, progeny, and the Licensed Materials including modified materials.

Between Genomeditech (Shanghai) Co., Ltd, and Licensee, Licensee is not permitted to modify cell lines in any way. The Licensee shall not share, distribute, sell, sublicense, or otherwise provide the Licensed Materials, or progenitors to third parties such as laboratories, departments, research institutions, hospitals, universities, or biotechnology companies for use other than for the purpose of outsourcing the Licensee's research.

Please refer to the Genomeditech Cell Line License Agreement for details.